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The Cary Eclipse fluorescence spectrophotometer was designed from the ground up, incorporating the latest optical and electronic technology. Included as standard are features such as corrected spectra, excitation and emission filters and extended range PMT detectors, so there won't be any nasty budget surprises.


Optional accessories such as the peltier thermostatted multicell holder and the microplate reader extend the capabilities of the instrument. With the flick of your wrist you can convert from using a standard cuvette to measuring a 384 well microplate.

Excellent sensitivity is a result of using an intense Xenon flash lamp, coupled with optimized grating blaze angles and coatings which ensure sensitivity across the whole wavelength range. Photosensitive samples are not exposed to continuous light as the Xenon flash lamp flashes only to acquire a data point.

Reduced volume cuvettes (40 μ L & 400 μ L) are available for use with the Cary Eclipse. The horizontal beam profile ensures excellent sensitivity even with such low volumes.

The large sample compartment makes it easy to install and remove accessories. You have plenty of room to fit your own research apparatus. We can even provide a baseplate for you to build upon.

The Schwarzschild collection optics capture a large portion of the light from the powerful Xenon flash lamp and direct it through the sample. This results in excellent sensitivity and low signal noise.

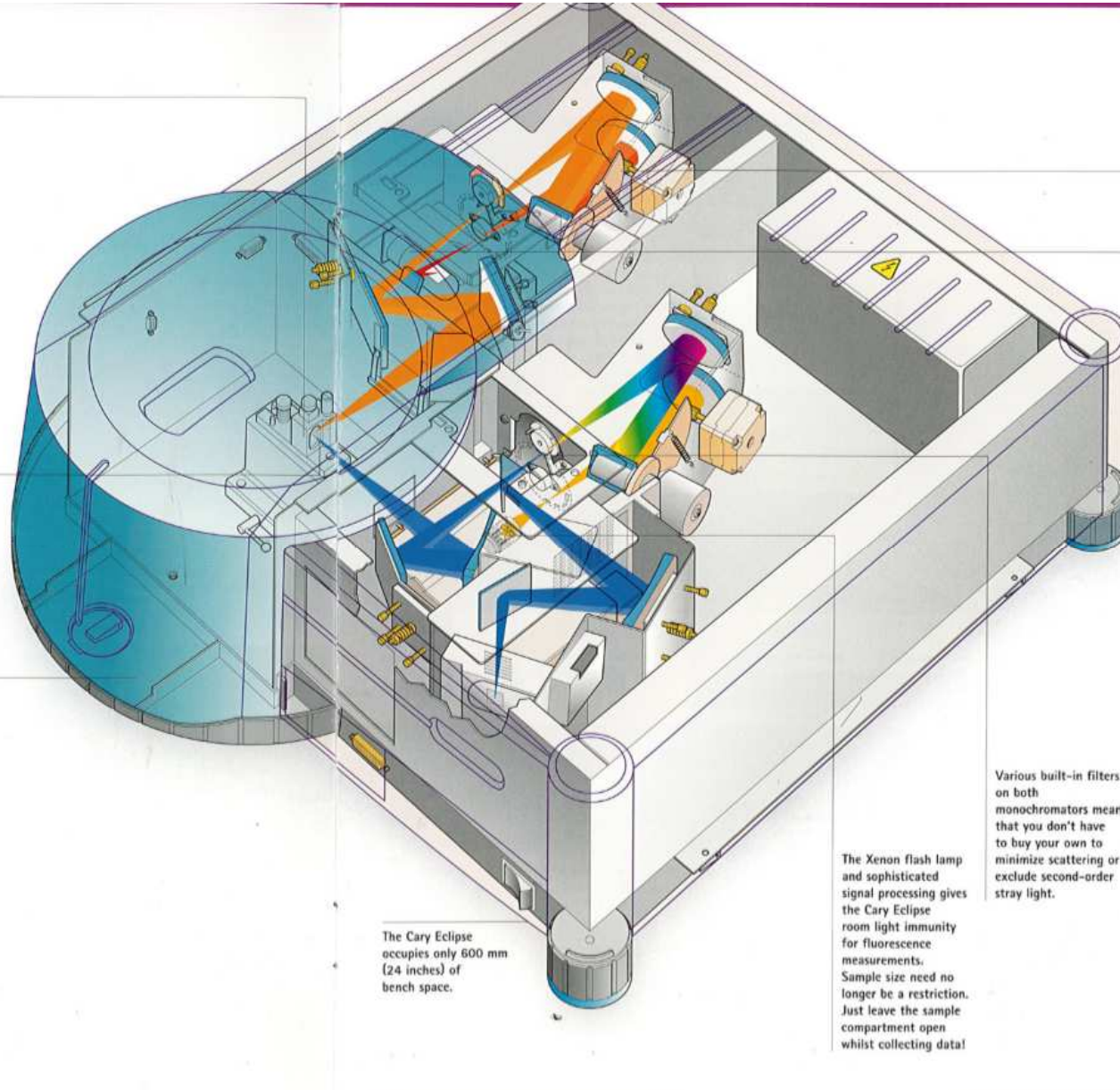Scan fast—the whole wavelength range in less than 3 seconds.

Red-sensitive Photomultiplier tube detectors extend sensitivity up to 900 nm without sacrificing UV performance.

Various built-in filters on both monochromators means that you don't have to buy your own to minimize scattering or exclude second-order stray light.

The Xenon flash lamp and sophisticated signal processing gives the Cary Eclipse room light immunity for fluorescence measurements. Sample size need no longer be a restriction. Just leave the sample compartment open whilst collecting data!

The Cary Eclipse occupies only 600 mm (24 inches) of bench space.